

# PDP Restart Procedure

To start PDP from scratch

1. Open NX session
  - Find NX icon at the top left of the screen. Click and select poltarg2p2.
  - The NX viewer is called "Remmina". If the icon is not in the right top corner, you can find it in the menu under internet.
2. Open Terminal
  - Once the NX window has finished opening, right click on the backdrop and select "terminal."
3. Start "System Start"
  - At terminal: "cd PDP"
  - "labview System\ Start.vi -- Configs/g2p.xml"
4. Start PDP
  - Click "Go!"
  - Wait until all windows open, and find "Display"
5. Select correct NMR channel and start taking data